**PROXECTO COMENIUS  
COMENIUS PROJECT**

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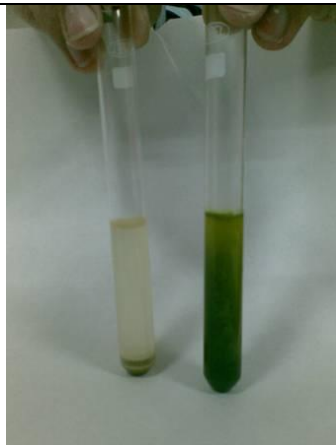
***Continuará...***

## Measurement of Iron concentration in galician plants.

### Procedure used to get the extracts

<p>1. Weigh a fixed amount of the vegetables and algae (18g in the first tests and 10g in the second one)</p>	
<p>2. Cut the vegetables in small pieces and we crush them by hand with the mixer or in the mortar (there are vegetables that are difficult to be handled with the electric mixer). The crushing operation is completed little by little adding about 25 cm<sup>3</sup> of water.</p>	
<p>3. Let them macerate until the following day, to be extract using a Büchner funnel (or using a vacuum pump). Separate the extract affluent from the filter paper.</p>	
<p>4. Prepare about 150 ml of tea infusion (using the traditional small bags). Two bags are enough and this tea infusion must rest during 24 hours before using.</p>	
<p>5. Distribute the extract in two probe tubes. Add the infusion in one and keep the other on as a pattern.</p>	
<p>6. For the vegetal and seaweed samples (approximately 10 ml of extract each one), add 6 ml of infusion tea with a pipette in each test tube.</p>	
<p>7. Let rest the solution and verifie the appearance of chemical precipitate in 30 minutes, 1 and 2 hours, and to the 24 hours.</p>	

## Observations Summary

Vegetable or seaweed	Observations 2 hours after adding “tea”	
<p><b>Espinaca (<i>Spinacia oleracea</i>)</b>  <b>Spinach</b>  <b>Epinard</b>  <b>(previous test)</b></p>	<p>Intense green precipitate (3 cm in the test tube)            No observed precipitate in the pattern sample</p>	
<p><b>Espárrago (<i>Asparagus officinalis</i>)</b>  <b>Asparagus</b>  <b>Asperge</b></p>	<p>Brown precipitate (0.5 cm).            No observed precipitate in the pattern sample</p>	
<p><b>Leituga (<i>Lactuca sativa</i>)</b>  <b>Lettuce</b>  <b>Laitue</b></p>	<p>Precipitate is not observed (turbid brown extract).            Pattern sample presents similar aspect.</p>	
<p><b>Grelos (<i>Brassica napus L.</i>)</b>  <b>Galician Turnip leaves</b>  <b>Feuilles de Navet (naveau, rave)</b></p>	<p>Intense green precipitate (4 cm in the test tube).            Green precipitate observed in the pattern sample also intense, but proportionally much smaller</p>	

<p><b>Repolo (<i>Brassica oleracea botrylis</i>)</b>  <b>Cabbage</b>  <b>Chou pomé</b></p>	<p>Intense green precipitate (3 cm in the tube).  Green precipitate observed in the pattern sample also intense, but proportionally much smaller</p>	
<p><b>Remolacha (<i>Beta vulgaris L.</i>)</b>  <b>Forage beet</b>  <b>Betteraves fourragères</b></p>	<p>Non observed precipitate (turbid green extract).  Pattern sample presents similar aspect.</p>	
<p><b>Verza (<i>Brassica oleracea botrylis</i>)</b>  <b>Cabbage leaves</b>  <b>Feuilles de Chou</b></p>	<p>Intense green precipitate (3,5 cm in the tube).  Green precipitate observed in the pattern sample also intense, but proportionally much smaller</p>	
<p><b>Canónigos (<i>Valerianella locusta</i>)</b>  <b>Lamb's salade</b>  <b>Mâche</b></p>	<p>Brown precipitate (2 cm).  Smaller Brown precipitate observed in the pattern sample</p>	

<p><b>Espinaca (<i>Spinacia oleracea</i>)</b> <b>Spinach</b> <b>Epinard</b></p>	<p>Intense green precipitate (3 cm in the test tube). No observed precipitate in the pattern sample.</p>	
<p><b>Leituga de mar (alga) (<i>Ulva rigida</i>)</b> <b>Sea lettuce (Seaweed)</b> <b>Laitue de mer (Algue)</b></p>	<p>Brown precipitate small particles (0.4 cm). No observed precipitate in the pattern sample</p>	

**Potassium hexacyanoferrate has been used to probe the presence of Fe(II) and Fe(III) in the samples.**

<b><i>Sample</i></b>	<b><i>Result immediately shown</i></b>
Ferrum (II) Chloride tetrahydrated (aq)	Blue intense
Ferrum (II) Amonium Sulfate hexahydrated (aq)	Blue intense
Ferrum (III) Nitrate (aq)	Brown (color tile)
Ferrum (III) Chloride tetrahydrated (aq)	Brown (tea color)
Spinach (extract)	No colour changed observed
Sea Lettuce (extract)	No colour changed observed

These results indicate that the concentration of ions are very small so we have to increase the concentration by heating the samples before repeating the tests.

Nevertheless, we have not carried out the test with the most of the samples, so maybe we get good results with the rest of plants and algae.

***To be continued...***